OPERATING INSTRUCTIONS

Humidity Cabinet (without Digital Data Recorder)

39-1600

| | | |
|---|---------------------|--|
| ELE International Chartmoor Road, Chartwell Business Park Leighton Buzzard, Bedfordshire, LU7 4WG England phone: +44 (0) 1525 249200 fax: +44 (0) 1525 249249 email: ele@eleint.co.uk http://www.ele.com ELE International, a division of Hach Lange Ltd. | Distributor: | ELE International Soiltest Product Division PO Box 389, Loveland, CO 80539 USA phone: +1 (800) 323 1242 fax: +1 (970) 663 9781 email: soiltest@eleusa.com http://www.eleusa.com |
| <p><i>In the interests of improving and updating its equipment, ELE reserves the right to alter specifications to equipment at any time.</i> ELE International 2018 ©</p> | | |

Contents

| | Section | Page |
|----------|--|-------------|
| 1 | General Description | 3 |
| 2 | Installation | 4 |
| 3 | Temperature Setting Procedure | 4 |
| 4 | Humidity Control | 7 |
| 5 | Maintenance | 7 |
| | APPENDIX A : Programming an Offset Parameter into a JUMO Dtron 316 Temperature Controller | 9 |
| | APPENDIX B : TLK38 Offset Instructions | 10 |
| | WEEE Directive | |

1 General Description

The cooled laboratory humidity cabinet provides the user with extremely accurate and safe temperature control over a wide temperature range.

1.1 Heating

Low wattage heaters are attached to the outer surface of the inner chamber in such a manner as to ensure even heating throughout the chamber.

1.2 Cooling

Direct expansion coils are coupled to a hermetic condensing unit via a solenoid valve. When the refrigeration switch is switched on, the compressor will run continuously. The coolant flow is controlled by a solenoid valve, which is electronically opened and closed by the temperature controller.

1.3 Temperature Controller

The temperature is accurately controlled by a Jumo dTRON 316 microprocessor temperature controller. The controller uses the signal from a PT100 sensor located in the chamber to control the current supplied to the heaters. The temperature can be easily adjusted. See Section 3 for details.

1.4 Over / Under Temperature Safety Cut Out

A TLK38 digital cut out protects the contents of your incubator by allowing you to program over & under temperature safety cut out set points. When the chamber temperature goes out of limits, an audible alarm and visible red neon will bring to your attention any problems. Reset buttons are provided below the cut out for you to manually reset the out-of-limit system.

1.5 Circulating Fan

Air is drawn up through the chamber by a circulating fan located in the top and returned to the base via a full width duct at the rear of the chamber. A door micro-switch stops the fan when the door is opened.

1.6 Humidity

A wide range of fixed relative humidities can be maintained by adding a suitable saturated salt solution to the base of the chamber. See Section 4 for details.

2 Installation

2.1 Connect the 3-core cable to a 240V AC, 50Hz mains supply as follows:

| | | |
|--------------|---|---------|
| BROWN | - | LIVE |
| BLUE | - | NEUTRAL |
| GREEN/YELLOW | - | EARTH |

2.2 The incubator is internally protected by a fuse, which isolates the cabinet from the electrical supply if excessive current is drawn. The fuse holder is located on the rear of the cabinet. The instruments on the control panel will illuminate when the supply is connected correctly and POWER switch is on.

2.3 General Precautions

A qualified electrician or other competent person must carry out any electrical work required to install the incubator.

Do not place the incubator in direct sunlight or near a heat source.

Make sure the incubator is not standing on its electrical supply cable.

Before the incubator is switched on for the first time, it should stand still for 2 hours after final positioning to allow the refrigerant to settle.

Keep all ventilation grilles clear. Failure to do so may result in component failure.

The refrigeration system gives off heat when operating. Always ensure there is adequate ventilation around the incubator.
NEVER PLACE ANYTHING ON TOP OF THE INCUBATOR.

Before any cleaning or maintenance work is carried out, the mains supply must be switched off and the plug removed from the electrical socket.

3 Temperature Setting Procedure

Switch both the POWER and COOLING switches ON. DO NOT switch the HIGH HUMIDITY switch on yet! *Please see Section 4.1 for more details.*

- 3.1 The JUMO dTRON 316 microprocessor temperature controller on your incubator has two LED displays. The upper display, which is red, represents the actual chamber temperature. The lower display, which is green, represents the target temperature. The controller has 4 buttons marked as follows;

- PGM** - Program button (Enters the programming mode).
- ▲** - Increases the temperature set point
- ▼** - Decreases the temperature set point.
- EXIT** - Exits the programming mode.

The **yellow 3** on the temperature controller shows when heat is being supplied. The **yellow 5** shows when cooling is being supplied.

- 3.2 The incubator operating temperature or set point (SP) can be altered by simply pressing either the **▲** button to increase the temperature or the **▼** button to decrease the temperature. When the required temperature has been entered, release all buttons and the green display will blink once to confirm that a new target temperature has been stored.

- 3.3 Over / Under Temperature Safety Cut Out

A TLK38 digital safety cut out unit provides essential over & under temperature protection for the products inside the chamber. The cut out has buttons labelled as follows:

- P** - Enters the programming mode.
- U** - *No function.*
- ▲** - **INCREASES** the displayed value.
- ▼** - **DECREASES** the displayed value.

3.4 Setting An Over Temperature Cut-Out Value

If, for example, you would like to operate your incubator at +20.0°C, an over temperature cut-out value needs to be programmed into the cut-out. We recommend that the cut-out value is set to **2 degrees higher** than your chamber temperature. In the example above, the over temperature cut-out set point needs to be set +22°C. To do this follow steps 1 - 3 below:

- a) Press the **[P]** button **once only**. Flashing **AL1** will be displayed plus a value (the **over temperature** cut out value).
- b) Press the **[▲]** button to increase the **over temperature** cut-out value or the **[▼]** button to decrease the **over temperature** cut-out value.
- c) Once the value of your choice is displayed, press the **[P]** button again and the new value will be stored and retained in permanent memory.

3.5 Over Temp Neon, Buzzer & Reset Button

If your incubator overheats, the buzzer will sound and the over temperature neon will illuminate. **The heaters will automatically disconnect to prevent further overheating.** Once the chamber temperature has stopped rising, and has fallen below the cut out set point, the heaters will re-connect automatically. The buzzer & neon will remain on until a member of staff has acknowledged the alarm. To reset the buzzer & alarm, press the reset button alongside the illuminated neon.

3.6 Setting An Under Temperature Cut Out Value

If, for example, you would like to operate your incubator at +20.0°C, an under temperature cut out value needs to be programmed into the cut out. We recommend that the cut-out value is set to **2 degrees lower** than your chamber temperature. In the example above, the under temperature cut out set point needs to be set +18°C. To do this follow steps 1 - 3 below:

- a) Press the **[P]** button **twice**. Flashing **AL2** will be displayed plus a value (the **under temperature** cut out value).
- b) Press the **[▲]** button to increase the **under temperature** cut out value or the **[▼]** button to decrease the **under temperature** cut out value.
- c) Once the value of your choice is displayed, press the **[P]** button again and the new value will be stored and retained in permanent memory.


3.7 Under Temp Neon, Buzzer & Reset Button

If your incubator overcools, the buzzer will sound and the under temperature neon will illuminate. **The cooling system will automatically disconnect to prevent further overcooling.** The buzzer & neon will remain on until a member of staff has acknowledged the alarm. To reset the buzzer & alarm, press the reset button alongside the illuminated neon.

- 3.8 Allow the chamber temperature to stabilise before putting your products into the chamber.

4 Humidity Control

A wide range of fixed and relative humidities can be achieved by adding a suitable saturated salt solution to the base of the chamber. The relative humidity of air in contact with saturated salt solutions reaches known equilibrium values at certain temperatures.

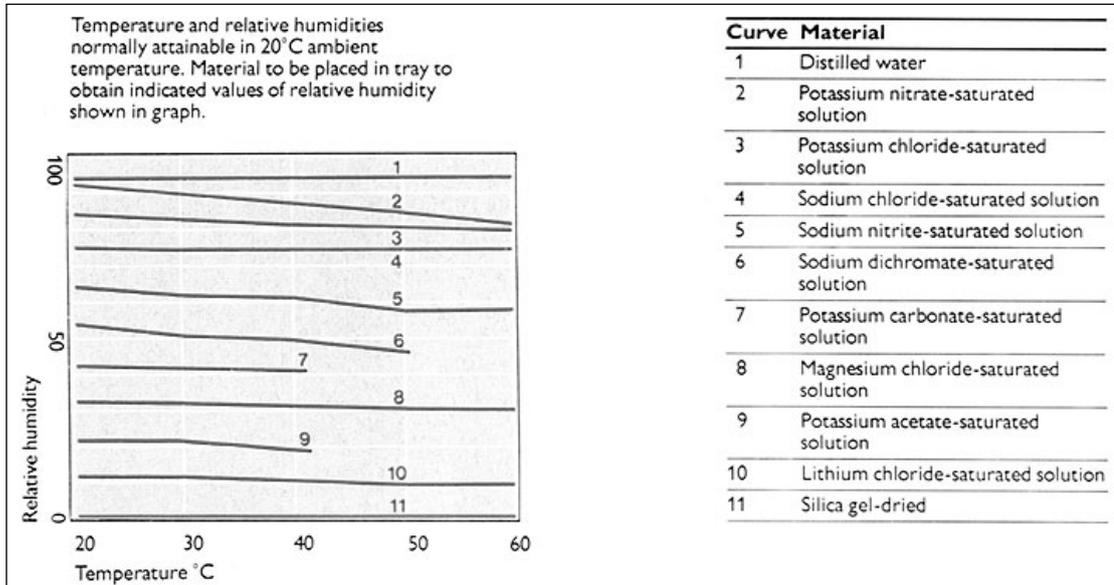Salts Graph

The graph above shows the temperatures and relative humidities for several of the more common saturated salts. The distilled water/saturated salt solution should be topped up regularly and should fully cover the floor AND high humidity element at all times.

4.1 High Humidity Option

A high humidity level of approx. 98%RH is maintained by placing distilled water in the base of the chamber and switching the HIGH HUMIDITY switch ON.

Do not use the HIGH HUMIDITY switch unless the floor of the chamber is fully covered with water. It is essential that the heating element on the chamber floor is covered at all times. This must be checked daily.

4.2 Allow the chamber humidity to stabilise before putting your products into the chamber.

5 Maintenance

No routine maintenance is necessary but regular cleaning of the inner chamber is essential.

5.1 Keep ventilation grilles clear and unobstructed.

- 5.2 The chamber should be regularly de-frosted as a substantial build-up of ice in the air circulation duct will restrict the airflow and affect the incubators performance. It is recommended that the service ports on the right of the incubator remain sealed with the threaded red bungs supplied to reduce the tendency for humidity in the air to condense out in the chamber.

APPENDIX A

Programming an Offset Parameter into a JUMO dTRON 316 Temperature Controller

If the temperature displayed by your replacement Jumo 316 controller does not match the actual chamber temperature, an offset parameter may have to be programmed into the controller. To do this, follow the steps below:

1. When the controller is in its normal mode, press the **PGM** button once.
2. **OPr** should be displayed.
3. Press the **▼** button twice. **ConF** should appear.
4. When **ConF** is displayed, press the **PGM** button once. **iNP** should appear.
5. Press the **PGM** button twice. **SEnS** should be displayed.
6. Press the **▼** button twice. **OFFS °C** should be displayed.
7. To adjust the offset value, press the **PGM** button (**OFFS** flashes), then use the **▲** button to increase the **OFFS** value, or the **▼** button to decrease the **OFFS** value. (If for example the temperature display was originally reading 2.0° too high, you need to put a negative offset value in of -2.0. If the temperature display was originally reading 1.0° too low, you need to put a positive offset value in of +1.0).
8. When the desired offset has been entered, **wait a few seconds** and the display will blink to confirm the new offset parameter has been stored.
9. Press the **EXIT** button four times.
10. The current chamber temperature will then be displayed which will include the additional offset value that you have just programmed in.
11. If at any stage during this procedure, the buttons are not pressed for 10 seconds or more, the controller will automatically return to its normal mode and you will have to restart from step 1.

APPENDIX B

TLK38 Offset Instructions

1. If the temperature displayed by the TLK38 does not match the calibrated test equipment, the TLK38 can be adjusted by means of an **offset parameter**.
2. Insert a calibrated probe (from your independent test equipment) into the incubator chamber via the access port.
3. Leave the incubator to reach its set point and stabilise.
4. Press and hold the **P** button for 5 seconds. **ConF** will be displayed.
5. Press the **P** button. **0** will be displayed.
6. Use the **▲** button to make the display read **38.1**
7. Press the **P** button. **5P** will be displayed.
8. Press the **▼** button once only. **1nP** will be displayed.
9. Press the **P** button. **5En5** will be displayed.
10. Press the **▼** button several times until **DFSt** is displayed.
11. Press the **P** button.
12. Change the value of **DFSt** from **0** (factory default) to whatever value is needed to correct the misreading display. (A positive or negative value can be entered).
13. Press the **P** button.
14. Hold the **▼** button for 1 second. **1nP** will be displayed.
15. Press the **▼** button. **DUt** will be displayed.
16. Press the **▼** button. **AL 1** will be displayed.
17. Press the **▼** button. **AL 2** will be displayed.
18. Press the **▼** button. **LbA** will be displayed.
19. Press the **▼** button. **rE9** will be displayed.
20. Hold the **▼** button for 2 seconds.
21. **The correct temperature should now be showing on the TLK38.**
22. POWER OFF / POWER ON to store the settings and reboot the incubator.

DO NOT CHANGE ANY OTHER SETTING IN THE TLK38 CONTROLLER

DIRECTIVE ON WASTE ELECTRICAL & ELECTRONIC EQUIPMENT (WEEE)

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